Arsenic (0–500 ppb) 2800000 (Arsenic Test Kit)

DOC326.97.00118

Test preparation

DANGER: ⚠ Hydrogen and arsine gasses are generated during the test. Work in a well-ventilated area away from open flames and other sources of ignition. Review Safety Data Sheets for safe handling, storage and disposal information.

CAUTION: ⚠ Review the Safety Data Sheets (MSDS/SDS) for the chemicals that are used. Use the recommended personal protective equipment.

- Do not expose reacted strips to direct sunlight. Reaction products are photosensitive and may turn dark.
- Do not allow test strips to touch the reaction vessel solution. Test strips react with gases, not solution.
- Orient the test strip pad **paper side down** and **centered** over the hole in the black cap so the generated gases can make good contact with the pad.
- Two samples may be analyzed simultaneously with this kit.
- This test measures inorganic arsenic. This test does not measure organic arsenic compounds, such as dimethylarsenic acid.
- Organic arsenic is a small proportion of the arsenic in most systems.
- To measure inorganic and organic arsenic (total arsenic), do the [Optional procedure to measure total arsenic](#) on page 1 before the test.

Interferences

Table 1 Interfering substances

Substance	Interference level
Sulfide	> 5 ppm
Selenium	> 1 ppm
Antimony	> 250 ppb
Tellurium	Likely to interfere, but not tested.

Table 2 Non-interfering substances

Substance	Highest concentration tested
Hardness	1000 ppm as CaCO ₃
Alkalinity	1000 ppm as CaCO ₃

Table 2 Non-interfering substances (continued)

Substance	Highest concentration tested
Iron	100 ppm
Temperature	10 to 40 °C (50 to 104 °F)

Required items


Description	Unit	Item no.
Arsenic Test Kit Reagent Set, includes: Arsenic test strips and arsenic Reagents #1, #2, #3, #4 and #5:	each	2799900
Arsenic test strips	100/pkg	—
Arsenic Reagent #1 Powder Pillows	100/pkg	—
Arsenic Reagent #2 Powder Pillows	100/pkg	—
Arsenic Reagent #3 Powder Pillows	100/pkg	—
Arsenic Reagent #4	250 g	—
Arsenic Reagent #5 Powder Pillows	100/pkg	—
Cap, Santoprene	2/pkg	4934800
Reaction vessel, 50-mL fill line	2/pkg	2800200
Scoop, 2 g	each	2799800

Optional procedure to measure total arsenic


Do the steps that follow before the test:

- Collect 50 mL of sample in a glass beaker.
- Add the contents of one Reagent #1 Powder Pillow to the sample. Swirl to dissolve.
- Add the contents of one Reagent #2 Powder Pillow to the sample. Swirl to dissolve. The solution may be cloudy.
- Put the beaker in a boiling water bath for 30 minutes.
- Put the prepared sample in the reaction bottle.
- Wait for the sample to decrease to room temperature.
- Do step 1 of the test procedure.
- Complete the test procedure. Start at step 6.

Test procedure—Arsenic (0, 10, 30, 50, 70, 300, 500 ppb)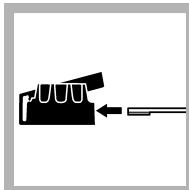

1. Insert a test strip into the cap so the pad completely covers the small opening. Close the flap and press to secure.

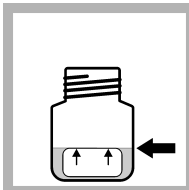
2. Fill the reaction bottle with sample to the fill line (50 mL).

3. Add one Reagent #1 Powder Pillow to the sample. Swirl slowly to dissolve.

4. Add one Reagent #2 Powder Pillow to the sample. Swirl slowly to dissolve. The solution may be cloudy.

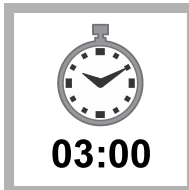
5. Wait 3 minutes (minimum).


6. Add one Reagent #3 Powder Pillow to the sample. Swirl slowly to dissolve. Some of the powder will not dissolve.


7. Wait 2 minutes (minimum). Swirl slowly to mix.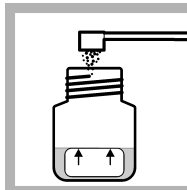


8. Use the plastic scoop to add 2 g of Reagent #4 to the sample. Swirl slowly to dissolve. Most of the powder will dissolve.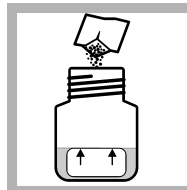


9. Add one Reagent #5 Powder Pillow to the sample.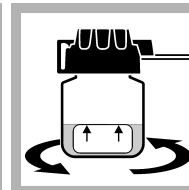

10. Immediately attach the cap to the reaction bottle. Swirl slowly for approximately 3 seconds. Do not shake or invert or allow sample to get on the strip.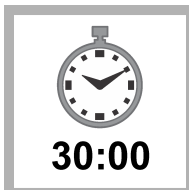

11. Wait 30 minutes, but no more than 35 minutes. Do not touch the reaction bottle for the full 30-minute reaction period.

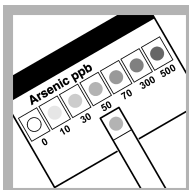12. Remove the test strip and immediately compare the developed color to the chart on the test strip bottle (0–500 ppb row). Read strips in the shade.

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