


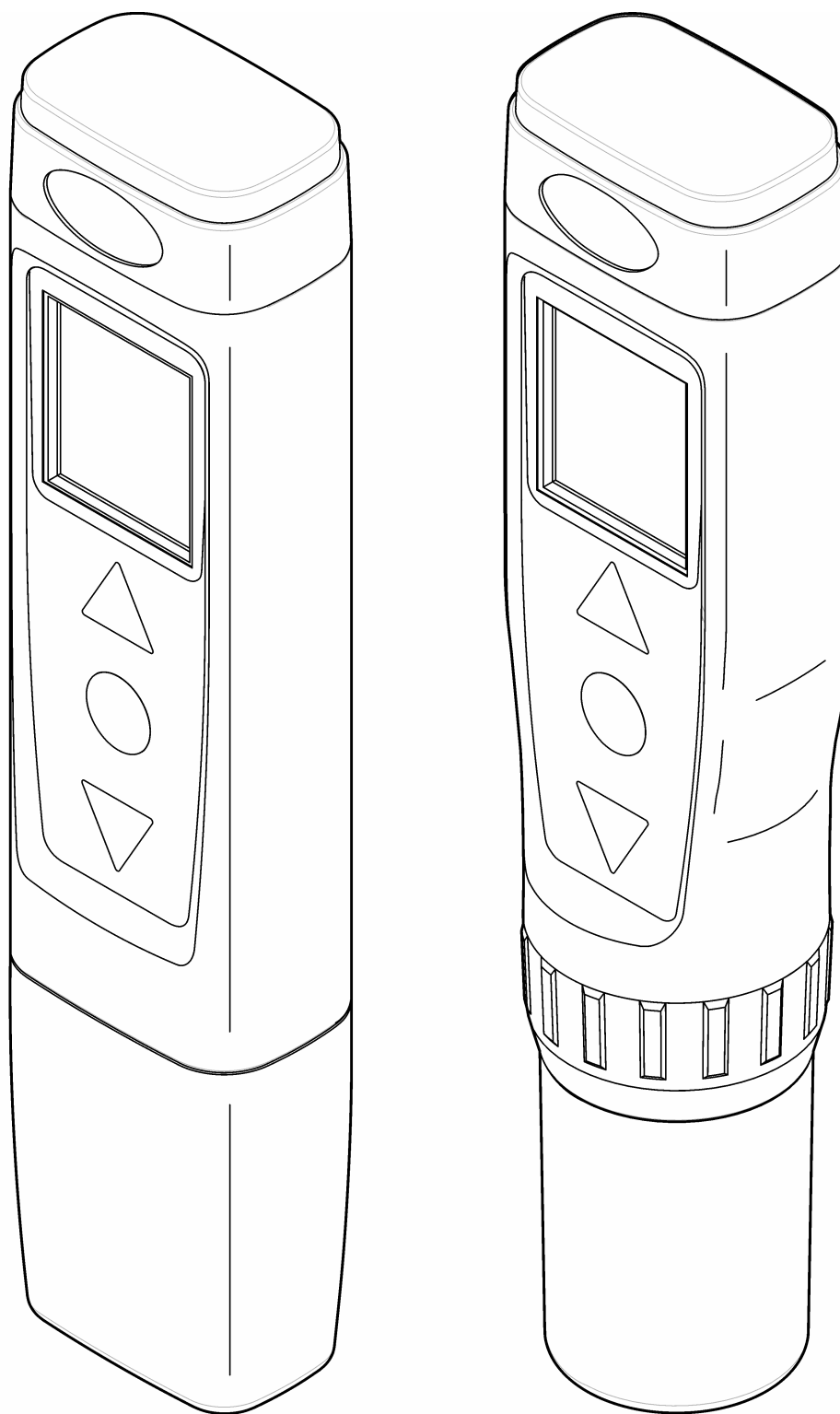

## Specifications

Specifications are subject to change without notice.

| Specification                          | Details                                                                                                                                                          |
|----------------------------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Dimensions (W x D x H)                 | 37 x 30 x 170 mm (1.45 x 1.18 x 6.69 in.)                                                                                                                        |
| Enclosure rating                       | IP67                                                                                                                                                             |
| Weight                                 | 135 g (0.297 lb) with batteries                                                                                                                                  |
| Power requirements (internal)          | AAA alkaline batteries (4x)                                                                                                                                      |
| Battery life                           | Pocket Pro: 450 hours<br>Pocket Pro+: 450 hours (200 hours with backlight on)                                                                                    |
| Operating temperature                  | 0 to 50 °C (32 to 122 °F)                                                                                                                                        |
| Storage temperature                    | –20 to 60 °C (–4 to 140 °F)                                                                                                                                      |
| Operating humidity                     | 80% (non-condensing)                                                                                                                                             |
| Altitude                               | 2000 m (6562 ft)                                                                                                                                                 |
| Instrument range                       | 0.0 to 14.0 pH                                                                                                                                                   |
| Resolution                             | Pocket Pro: 0.1 pH; Pocket Pro+: 0.01 pH                                                                                                                         |
| Accuracy <sup>1</sup>                  | Pocket Pro: ±0.1 pH; Pocket Pro+: ±0.02 pH                                                                                                                       |
| Calibration points                     | Pocket Pro: 3 points (auto); Pocket Pro+: 3 points (auto) or 2 points (custom)<br><b>Note:</b> Temperature calibration is not available.                         |
| Auto-recognition calibration standards | USA: 4.01, 7.00, 10.01 pH; NIST: 4.01, 6.86, 9.18 pH                                                                                                             |
| Certifications                         | CE mark, FCC, Industry Canada, KC Mark, RCM, China RoHS                                                                                                          |
| Warranty                               | 6 months for Pocket Pro tester, 1 year for Pocket Pro+ tester and 6 months for replacement sensor for manufacturing faults only. Damage from use is not covered. |

## General information

In no event will the manufacturer be liable for direct, indirect, special, incidental or consequential damages resulting from any defect or omission in this manual unless otherwise required by applicable law or contract between the parties. The manufacturer reserves the right to make changes in this manual and the products it describes at any time, without notice or obligation. Revised editions are found on the manufacturer's website.

## Safety information

The manufacturer is not responsible for any damages due to misapplication or misuse of this product including, without limitation, direct, incidental and consequential damages, and disclaims such damages to the full extent permitted under applicable law. The user is solely responsible to identify critical application risks and install appropriate mechanisms to protect processes during a possible equipment malfunction.

Please read this entire manual before unpacking, setting up or operating this equipment. Pay attention to all danger and caution statements. Failure to do so could result in serious injury to the operator or damage to the equipment.

If the equipment is used in a manner that is not specified by the manufacturer, the protection provided by the equipment may be impaired. Do not use or install this equipment in any manner other than that specified in this manual.

---

<sup>1</sup> Based on a 3-point calibration and the calibration standards at the same temperature as the samples measured. Also valid for 5.5 to 8.5 pH based on a 1-point calibration, 0.0 to 8.5 pH based on a 2-point calibration with pH 7 and pH 4 standards, or 5.5 to 14 pH based on a 2-point calibration with pH 7 and pH 10 standards.

## Use of hazard information

### DANGER

Indicates a potentially or imminently hazardous situation which, if not avoided, will result in death or serious injury.

### WARNING

Indicates a potentially or imminently hazardous situation which, if not avoided, could result in death or serious injury.

### CAUTION

Indicates a potentially hazardous situation that may result in minor or moderate injury.

### NOTICE

Indicates a situation which, if not avoided, may cause damage to the instrument. Information that requires special emphasis.

## Precautionary labels

Read all labels and tags attached to the instrument. Personal injury or damage to the instrument could occur if not observed. A symbol on the instrument is referenced in the manual with a precautionary statement.



This symbol, if noted on the instrument, references the instruction manual for operation and/or safety information.



Electrical equipment marked with this symbol may not be disposed of in European domestic or public disposal systems. Return old or end-of-life equipment to the manufacturer for disposal at no charge to the user.

## Certification

### Canadian Radio Interference-Causing Equipment Regulation, ICES-003, Class B:

Supporting test records reside with the manufacturer.

This Class B digital apparatus meets all requirements of the Canadian Interference-Causing Equipment Regulations.

Cet appareil numérique de classe B répond à toutes les exigences de la réglementation canadienne sur les équipements provoquant des interférences.

### FCC Part 15, Class "B" Limits

Supporting test records reside with the manufacturer. The device complies with Part 15 of the FCC Rules. Operation is subject to the following conditions:

1. The equipment may not cause harmful interference.
2. The equipment must accept any interference received, including interference that may cause undesired operation.


Changes or modifications to this equipment not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment. This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference, in which case the user will be required to correct the interference at their expense. The following techniques can be used to reduce interference problems:

1. Move the equipment away from the device receiving the interference.
2. Reposition the receiving antenna for the device receiving the interference.
3. Try combinations of the above.

Product overview

The Pocket Pro™ pH tester and Pocket Pro™+ pH tester measure the pH of general water samples. Refer to [Figure 1](#). These testers are waterproof and float.


Figure 1 Product features

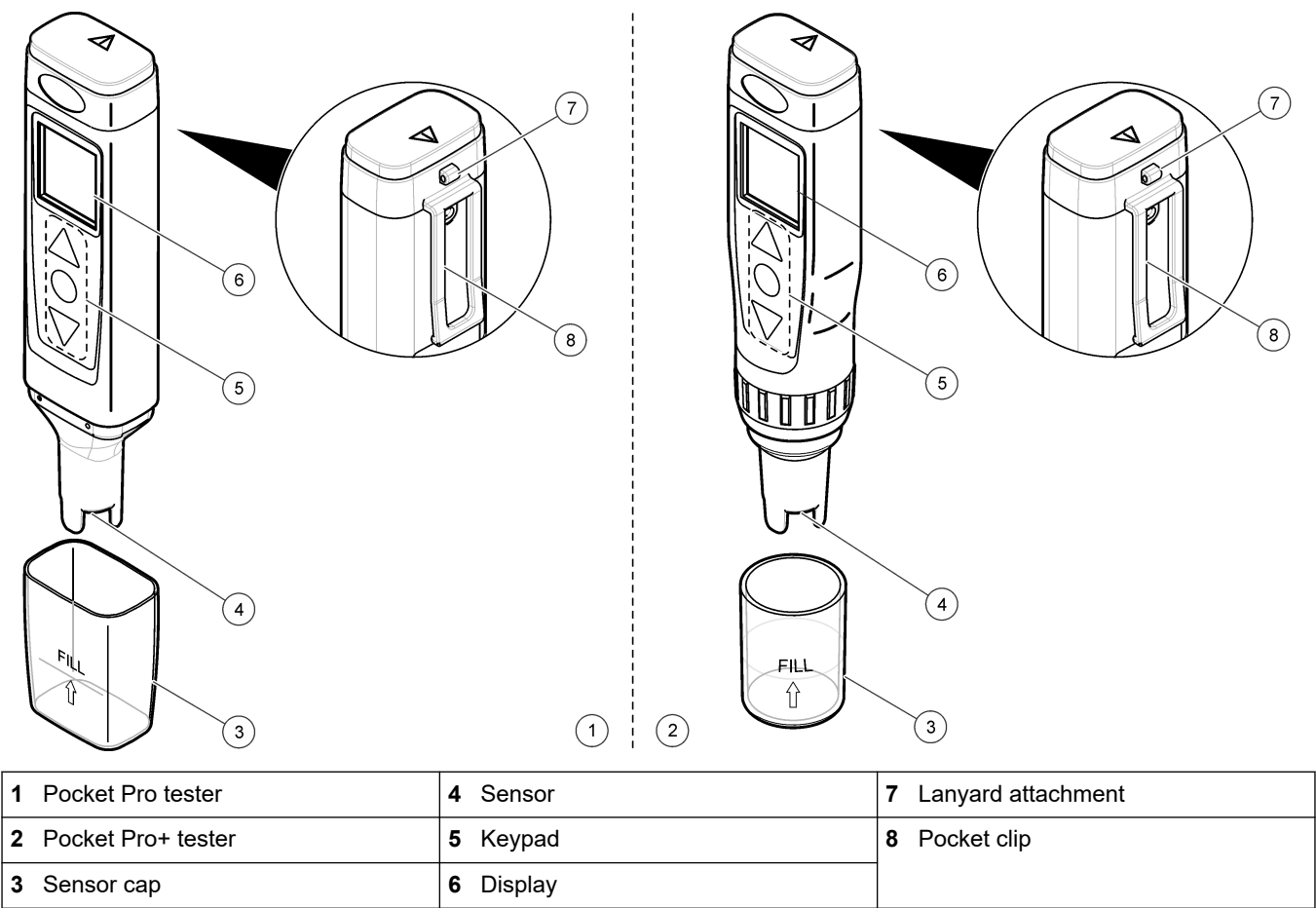

Product components

Make sure that all components have been received. Refer to [Figure 2](#). If any items are missing or damaged, contact the manufacturer or a sales representative immediately.

Figure 2 Product components

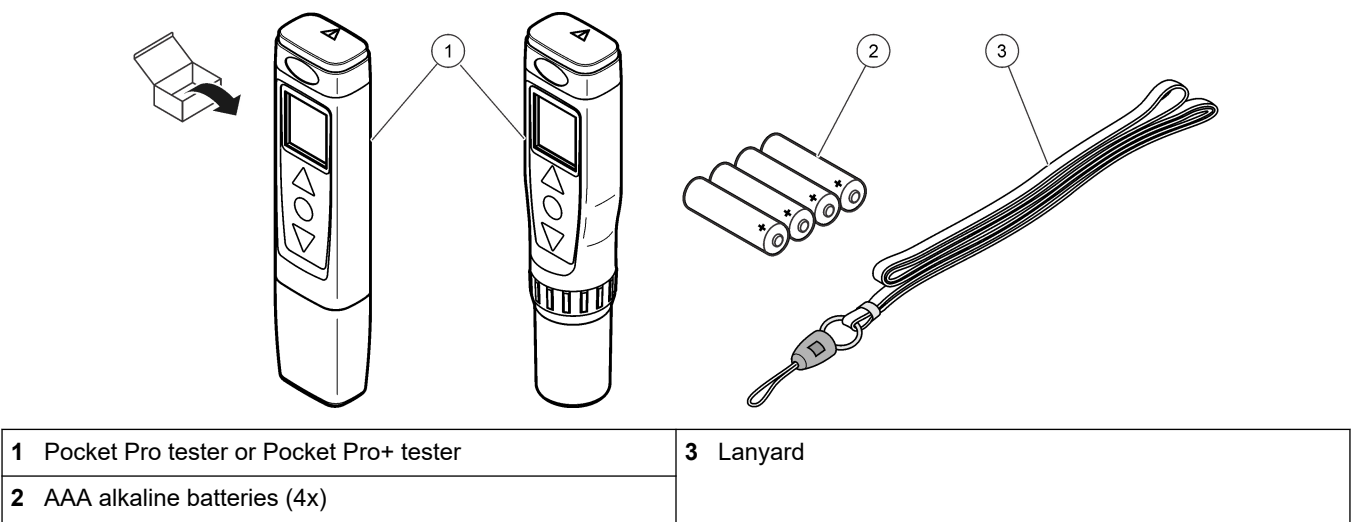

## Install the batteries

### ⚠ CAUTION



Explosion hazard. Incorrect battery installation can cause the release of explosive gases. Be sure that the batteries are of the same approved chemical type and are inserted in the correct orientation. Do not mix new and used batteries.


### ⚠ WARNING



Fire hazard. Battery substitution is not permitted. Use only alkaline batteries.

Install the four AAA alkaline batteries in the tester. Refer to the illustrated steps in [Figure 3](#).

**Figure 3** Install the batteries

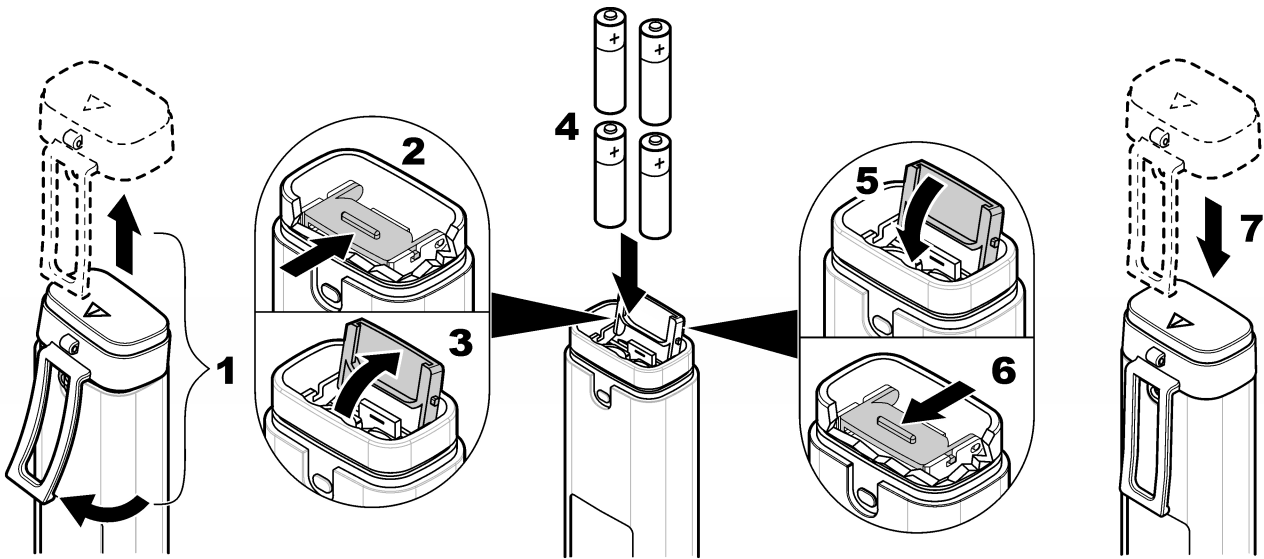

# User interface and navigation

## Display description

Figure 4 shows the measurements, calibration standard information and indicator icons shown on the display. Table 1 gives descriptions of the icons.

Figure 4 Display overview

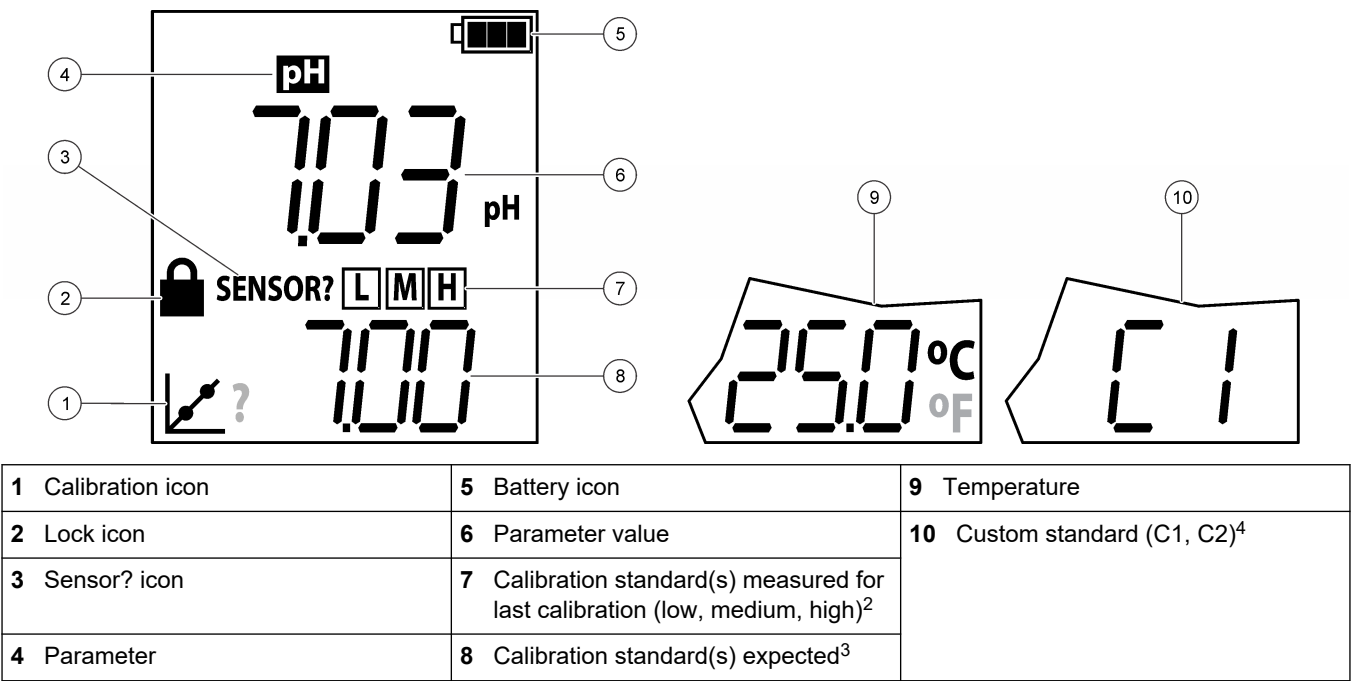


Table 1 Display icons


| Icon    | Description                                                                                                                                                                                                                                                                     |
|---------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|         | Shows the battery power level. Flashes when the battery power is less than 10%.                                                                                                                                                                                                 |
|         | Shows when the lock feature is on. When the lock feature is on, the parameter value on the display does not change. Push  to set the lock feature to on or off.                                                                                                                 |
| Sensor? | Refer to <a href="#">Troubleshooting</a> on page 10.                                                                                                                                                                                                                            |
|         | Shows when the tester is in calibration mode or when a calibration is due. If "?" shows next to the calibration icon, the last calibration was not successful. When the ACAL setting is set to Yes (default), "?" shows next to the calibration icon when a calibration is due. |

## Keypad description

Figure 5 and Figure 6 show the Pocket Pro and Pocket Pro+ keypads. Table 2 gives the key descriptions.


<sup>2</sup> Shows after auto calibration  
<sup>3</sup> Shows during auto calibration  
<sup>4</sup> Shows during custom calibration.

**Figure 5 Keypad overview - Pocket Pro**

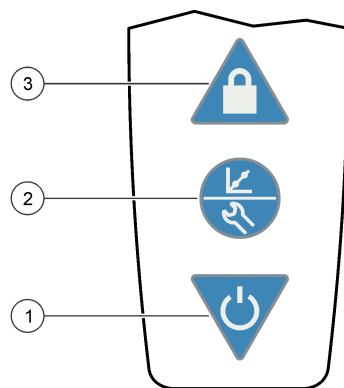

|             |                            |            |
|-------------|----------------------------|------------|
| 1 Power key | 2 Calibration/Settings key | 3 Lock key |
|-------------|----------------------------|------------|

**Figure 6 Keypad overview - Pocket Pro+**

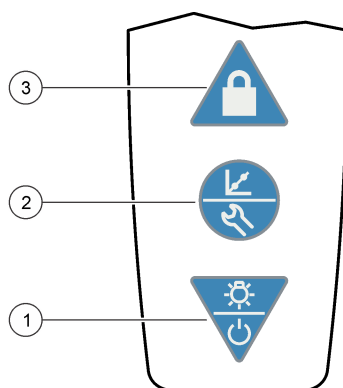

|                       |                            |            |
|-----------------------|----------------------------|------------|
| 1 Power/Backlight key | 2 Calibration/Settings key | 3 Lock key |
|-----------------------|----------------------------|------------|

**Table 2 Key functions**

| Key | Description                                                                                                                                                                                                                                                                                                                                                                   |
|-----|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|     | Push and hold to set the power to on or off.                                                                                                                                                                                                                                                                                                                                  |
|     | Push and hold to set the power to on or off.<br>Push to set the backlight to on or off. After 1 minute of no activity, the backlight switches off.                                                                                                                                                                                                                            |
|     | Push to start a calibration. To exit a calibration, push and hold.<br>Push and hold until "SEt" shows on the display to go to the settings menu. To exit the settings menu, push and hold until "End" shows on the display. When in the settings menu, push to scroll through the settings.<br><b>Note:</b> Power cannot be set to off while in settings or calibration mode. |
|     | Push to set the lock feature to on or off. When the lock feature is on, the lock icon shows and the parameter value on the display does not change.                                                                                                                                                                                                                           |

## Calibration

### ⚠ CAUTION



Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.

Calibrate the tester before initial use and when:

- Results drift

- 
- Results are not accurate
  - "?" shows next to the calibration icon

Calibrate the tester with the:

- Auto calibration procedure
- Custom calibration procedure<sup>5</sup>

Before the initial calibration and after a dry storage, soak the sensor for several minutes in the sample or tap water.

## Auto calibration

**Items to collect:** One, two or three auto-recognition calibration standards

1. Set the power to on.
2. Remove the cap from the sensor.
3. Push  to go to calibration mode.  
The auto-recognition standard(s) to measure shows on the bottom line.  
*Note: If "C1" shows on the bottom line, do not continue. Set the tester to auto calibration mode. Refer to [Configure the settings](#) on page 9.*
4. Rinse the sensor and cap with deionized water and blot dry.
5. Pour the auto-recognition standard shown into the cap to the fill line.
6. Put the sensor fully into the cap.
7. When the measurement is stable, push  to save the measurement.  
The measured value flashes three times.
8. Optional: To measure another calibration standard (maximum of 3), do steps 4–7 again.
9. Push and hold  to go to continuous measurement mode. "END" shows on the display.  
*Note: "ECAL" shows on the display if the calibration was not successful. Refer to [Troubleshooting](#) on page 10.*
10. Rinse the sensor and cap with deionized water and blot dry.

## Custom calibration (Pocket Pro+ only)

**Items to collect:** One or two calibration standards or samples of known pH value

1. Set the power to on.
2. Remove the cap from the sensor.
3. Go to the settings menu. Make sure that the bUFR setting is set to Cus (custom standard). Refer to [Configure the settings](#) on page 9.
4. Push  to go to calibration mode.  
"C1" shows on the bottom line.
5. Rinse the sensor and cap with deionized water and blot dry.
6. Pour the calibration standard or sample into the cap to the fill line.
7. Put the sensor fully into the cap.
8. When the measurement is stable, push  and  until the pH value of the calibration solution or sample shows on the display.
9. Optional: To measure a second pH calibration standard or sample of known value, push , then do steps 5–8 again.
10. When the last measurement is stable, push and hold  to save the calibration and go to continuous measurement mode. "END" shows on the display.  
*Note: "ECAL" shows on the display if the calibration was not successful. Refer to [Troubleshooting](#) on page 10.*
11. Rinse the sensor and cap with deionized water and blot dry.

---

<sup>5</sup> Custom calibration can only be done on the Pocket Pro+ tester.

## Measurement

### NOTICE

For storage, do not keep the sensor in a solution or damage to the sensor can occur.

**Note:** Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. Shake the tester from side to side to remove air bubbles.

1. Set the power to on.
  2. Remove the cap from the sensor.
  3. Before the measurement and after a dry storage, soak the sensor for several minutes in the sample or tap water.
  4. If the lock icon shows on the display, push  to go to continuous measurement mode.
  5. Rinse the sensor and cap with deionized water and blot dry.
  6. Pour the water sample into the cap to the fill line.
  7. Put the sensor fully into the cap. The measured value shows on the top line.
  8. To keep the measured value on the display when the sensor is removed from the sample, push .
- Note:** The lock icon shows on the display when the measurement is stable.
9. To measure another sample, do steps 4–8.
  10. When done with measurements:
    - a. Rinse the sensor and cap with deionized water.
    - b. Put the cap on the tester.
    - c. Set the power to off.
    - d. (Optional) Add some drops of tap water or deionized water to the storage cap to keep the sensor prepared for use.
    - e. Install the storage cap on the sensor.

## Advanced operation

### Configure the settings

1. Push and hold  until "SEt" shows on the display.
2. Push  to scroll through the settings. The current value of the setting shows on the bottom line.

| Option      | Description                                                                                                                                                                                                        |
|-------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>Unit</b> | Select the temperature unit that shows on the display— C (Celsius) or F (Fahrenheit).                                                                                                                              |
| <b>bUFr</b> | Select the pH buffers that are used for auto calibration—USA (4.01, 7.01, 10.01, default), NIST (4.01, 6.86, 9.18) or Cus (custom standard).<br><b>Note:</b> "Cus" can only be selected on the Pocket Pro+ tester. |
| <b>AOFF</b> | Set the auto-off feature to On (enable, default) or Off (disabled). When set to On, power is automatically set to off after 8 minutes of no activity.                                                              |
| <b>ACAL</b> | Enable or disable pH calibration reminder—Yes (enable, default) or No (disable). When set to Yes, "?" shows next to the calibration icon on the pH screen when a pH calibration is due.                            |
| <b>rSEt</b> | Change the settings to the factory defaults—Yes or No (default). When set to Yes, changes the settings to the factory settings and default values.                                                                 |

3. To change the value of the setting, push  or .
4. When done with changes, push and hold  until "End" shows to go to continuous measurement mode.

## Maintenance

### ⚠ CAUTION



Multiple hazards. Only qualified personnel must conduct the tasks described in this section of the document.

### Clean the sensor

Clean the sensor when:

- "SENSOR?" shows on the display
- Stabilization is slow
- Results drift or are not accurate
- Calibration failure occurs

1. Soak the sensor in the applicable cleaning agent. Refer to [Table 3](#).
2. Rinse or soak the sensor in deionized water for 1 minute.

**Table 3 Cleaning agents**

| Contaminant           | Cleaning agent                       | Time              |
|-----------------------|--------------------------------------|-------------------|
| Grease, oils and fats | Electrode cleaning solution          | 2 hours maximum   |
| Mineral buildup       | 10% hydrochloric acid (HCl) solution | 5 minutes maximum |

### Replace the batteries

When the battery icon flashes or the tester will not come on, replace all four batteries. Refer to [Install the batteries](#) on page 5.

### Replace the sensor

**Note:** Only Pocket Pro+ testers have replaceable sensors.

To replace the sensor, refer to the instructions supplied with the sensor.

## Troubleshooting

| Message                    | Possible cause                                                                                                                                                                                                     | Solution                                                                                                                                                                                                                                                                              |
|----------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| SENSOR?                    | The calibration slope is $\pm 10$ –15%.                                                                                                                                                                            | Gently clean the sensor. Refer to <a href="#">Clean the sensor</a> on page 10.                                                                                                                                                                                                        |
| ECAL                       | Calibration failure. The pH calibration slope is greater than $\pm 15$ %.                                                                                                                                          | Gently clean the sensor. Refer to <a href="#">Clean the sensor</a> on page 10. Calibrate again. If calibration failure continues, replace the tester or sensor as applicable.                                                                                                         |
| "- - - -"<br>(top line)    | The parameter value is not within the measurement range of the tester. Refer to <a href="#">Specifications</a> on page 2.                                                                                          | Make sure that the value of the sample is within the measurement range of the tester. Gently clean the sensor. Refer to <a href="#">Clean the sensor</a> on page 10. Calibrate as necessary.                                                                                          |
| "- - - -"<br>(bottom line) | The temperature value is not within the operating temperature range of the tester or a temperature sensor failure occurred. Refer to <a href="#">Specifications</a> on page 2 for the operating temperature range. | Make sure that the sample temperature is within the operating temperature range of the tester. Contact technical support as necessary.<br><br><b>Note:</b> The tester can still be used if a temperature sensor failure has occurred, but without automatic temperature compensation. |
| Battery icon flashes       | The batteries have less than 10% power remaining.                                                                                                                                                                  | Replace all four batteries. Refer to <a href="#">Install the batteries</a> on page 5.                                                                                                                                                                                                 |

## Replacement parts and accessories

### WARNING



Personal injury hazard. Use of non-approved parts may cause personal injury, damage to the instrument or equipment malfunction. The replacement parts in this section are approved by the manufacturer.

**Note:** Product and Article numbers may vary for some selling regions. Contact the appropriate distributor or refer to the company website for contact information.

**Table 4 Replacement parts**

| Description            | Quantity | Item no. |
|------------------------|----------|----------|
| AAA alkaline batteries | 4/pkg    | 4674300  |
| Lanyard                | 1        | 201305   |
| SINGLET, 7.00 pH       | 20/pkg   | 2770120  |
| pH sensor, replacement | 1        | 9532001  |

**Table 5 Accessories**

| Description                 | Quantity | Item no. |
|-----------------------------|----------|----------|
| SINGLET, 4.01 pH            | 20/pkg   | 2770020  |
| SINGLET, 10.01 pH           | 20/pkg   | 2770220  |
| Electrode cleaning solution | 500 mL   | 2965249  |
| Hydrochloric Acid (HCl)     | 2.5 L    | 13406    |
| Hydrochloric Acid (HCl)     | 500 mL   | 13449    |

---

**HACH COMPANY World Headquarters**

P.O. Box 389, Loveland, CO 80539-0389 U.S.A.  
Tel. (970) 669-3050  
(800) 227-4224 (U.S.A. only)  
Fax (970) 669-2932  
orders@hach.com  
www.hach.com

**HACH LANGE GMBH**

Willstätterstraße 11  
D-40549 Düsseldorf, Germany  
Tel. +49 (0) 2 11 52 88-320  
Fax +49 (0) 2 11 52 88-210  
info-de@hach.com  
www.de.hach.com

**HACH LANGE Sàrl**

6, route de Compois  
1222 Vérenaz  
SWITZERLAND  
Tel. +41 22 594 6400  
Fax +41 22 594 6499

