

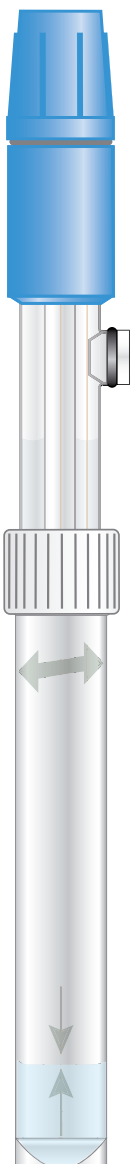
52 21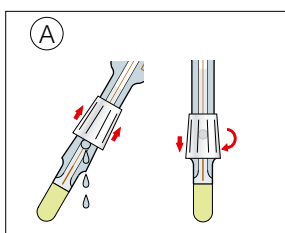

DOC022.98.90542
09-2015, Edition 1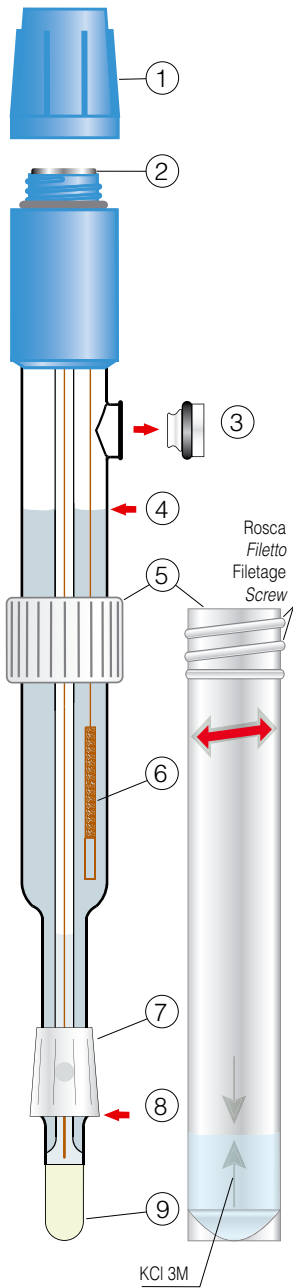

Puntos clave. *Punti chiave.* Points clé. *Key points.*

- ① Tapón protector.
Tappo protettore.
Bouchon protecteur.
Protective cap.
- ② Cabezal con conector S7.
Testa con connettore S7.
Tête avec connecteur S7.
S7 connector head.
- ③ Orificio de relleno de electrolito.
Orificio di riempimento per l'elettrolita.
Orifice de remplissage de l'électrolyte.
Electrolyte filling port.
- ④ Electrolito de referencia.
Elettrolita di riferimento.
Électrolyte de référence.
Reference electrolyte.
- ⑤ Protector de almacenamiento.
Protettore per la conservazione.
Protecteur de stockage.
Storage protector.
- ⑥ Elemento de referencia.
Elemento di riferimento.
Élément de référence.
Reference element.
- ⑦ Diafragma esmerilado.
Diaframma smerigliato.
Diaphragme à rodage.
Sleeve diaphragm.
- ⑧ Nivel mínimo de muestra.
Livello minimo di campione.
Niveau minimum d'échantillon.
Minimum sample level.
- ⑨ Membrana sensible.
Membrana sensibile.
Membrane sensible.
Sensitive glass membrane.

Electrodo de pH 52 21

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pH electrode 52 21

Introduction

The **52 21** is a pH electrode with S7 screw cap, suitable for difficult samples.

Electrode preparation

Unscrew the storage protector ⑤.

Verify that there are not air bubbles inside the membrane. They can be removed by shaking the electrode as a clinical thermometer.

Remove the cap from the electrolyte filling port ③. Before measuring, slide the plastic ring and return it to its original position, by pressing softly, to ensure the electrical contact, see fig A. If it is necessary, refill with reference electrolyte.

Calibration

⚠ See *pH-meter instruction manual*.

Measurements. Recommendations

- During calibration and measurement the buffers and the samples must be slightly stirred. Identical stirring speed in both cases.
- The sample must cover the electrode's diaphragm ⑦.
- After measurement, immediately rinse the electrode with distilled water. Dry it with soft tissue, only by contact, without rubbing.
- Between measurements the electrode's membrane should be immersed in aqueous solution. The most appropriate place is the storage protector with some electrolyte.

⚠ *Do not leave immersed in distilled water.*

Maintenance

- Periodically refill the electrode with electrolyte. Maintain the level as high as possible ④.
- Maintain the membrane's surface clean ⑨. Protect it from friction or knocking.
- When an electrode gets contaminated, the simple rinsing with distilled water is not sufficient. The most suitable cleaning solution is one which acts in a more selective way over the contamination (look for HACH regeneration solutions).
- **Re-hydration.** A "dry" electrode recovers by immersing it in HACH cleaning electrode solution or in diluted HCl solution for several hours. It is recommended to read the information supplied with the electrode's regenerating solutions.

Troubleshooting

... the meter always measures the same value.

Short circuit in the electrode or connector.

Substitute the electrode.

... unstable reading.

Air bubble in the membrane.

Shake the electrode as a thermometer.

... impossible to perform calibration.

... very slow response.

Dirt on the diaphragm or the membrane.

De-hydrated membrane.

Electrode ageing or wear.

Clean, regenerate, substitute electrode.

Storage

The connector should be protected from humidity by the protective cap ①.

Place the electrode in the storage protector ⑤ with KCl 3M.

Close the refilling port with the stopper ③.

Life of an electrode

The average life expectancy of a pH electrode is approximately one year depending on the working conditions.

An electrode used for more than six months that does not response appropriately after regeneration, must be replaced by a new one.

To avoid accidental breakage of the glass body electrodes, the use of the working protector, code 91 62, is recommended.

 *If an electrode has too short lifespan, less than three months, ask us about it. Probably there is a more suitable electrode for the application.*

Specifications

Measuring range	0 ... 12 pH
Operating temp.	0 ... 60 °C
Reference element	encapsulated Ag/AgCl crystals
Diaphragms	sleeve
Electrolyte	KCl 3M
Body material	glass
Min. immersion depth	20 mm 

Applications

- Samples with very low conductivity.
- Distilled water.
- Samples with colloids, wines, paints, etc.
- Very viscous sample, emulsions, creams...

Limitations

Temperatures > 60 °C.

Warranty

HACH guarantees this electrode against manufacturing defects.

Validity: 6 months from shipment date.

Limitations:

- Accidental damage.
- Inadequate applications.
- Non-fulfilment of the recommendations.
- Normal wear and tear of the electrode.

QUALITY CERTIFICATE

The attached electrode **52 21** has passed the quality test and follows the next specifications:

- Asymmetry potential < ± 15 mV.
- Sensitivity, pH 4...7 (at 25 °C) > 98%.
- Response time, pH 4...7 < 20 s.

HACH COMPANY World Headquarters

P.O. Box 389, Loveland, CO 80539-0389 U.S.A.
Tel. (970) 669-3050
(800) 227-4224 (U.S.A. only)
Fax (970) 669-2932
orders@hach.com
www.hach.com

HACH LANGE GMBH

Willstätterstraße 11
D-40549 Düsseldorf, Germany
Tel. +49 (0) 2 11 52 88-320
Fax +49 (0) 2 11 52 88-210
info@de.hach.com
www.de.hach.com

HACH LANGE Sàrl

6, route de Compois
1222 Vérenaz
SWITZERLAND
Tel. +41 22 594 6400
Fax +41 22 594 6499